

Air compressor filter kit June 2025


The SRI air compressor replacement kit (8670-1070) comes with a particle filter and 16 inches of 1/4" x 3/8" PVC tubing to connect the air compressor outlet to the filter. Also included are 3 crimp connectors to connect the white and black power wires.

See the you-tube video about how to replace the built-in air compressor on the SRI you-tube channel [youtube.com@SRIGCTraining/](https://www.youtube.com/@SRIGCTraining/) playlists

Most SRI GCs will already be equipped with a filter (making it easy to replace) but if there is no existing filter the downstream side of the filter will need to be connected to the existing tubing pathway with some smaller PVC tubing included in the kit.

The original air compressor may have a metal tube connected to its output

The replacement air compressor has a barb fitting which the PVC slips over and is secured by a clamp.

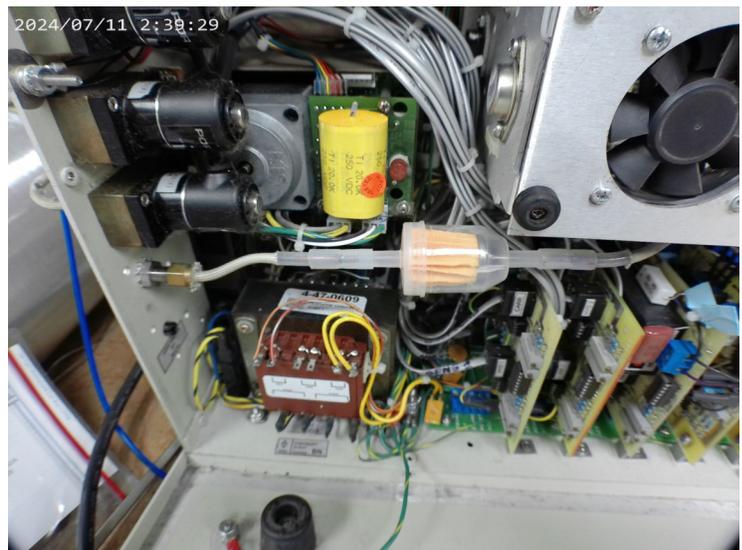

SRI Tech Support:
310-214-5092