


Moving and Packing the SRI 8610C,D or V and 310C GC for shipping

ENSURE that all 6 screws holding the bottom cover are fastened.

Lift the 8610C&D model GCs as shown to right.

Lift the Model 310C GC as shown.


Lift the Model 8610V as shown

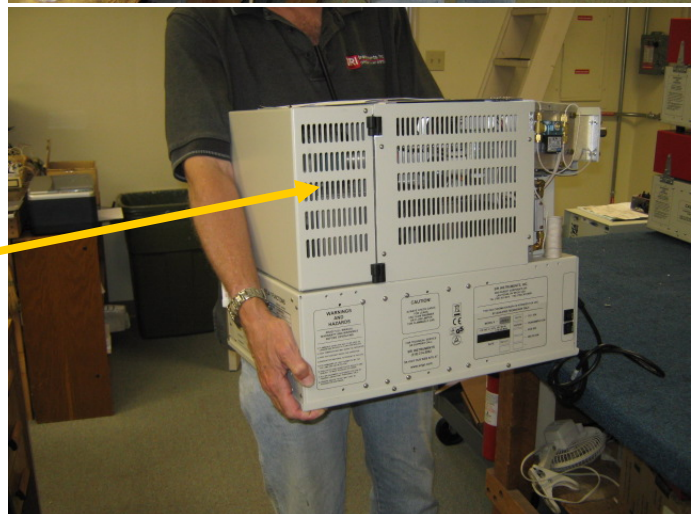

Moving and Packing the SRI 8610C,D or V and 310C GC for shipping

The photos in this document show the 8610C GC, but the 310C GC is very similar except a little smaller in size.

Place the SRI 8610C GC on the blue foam. The foam has holes for the GC feet to sit in.

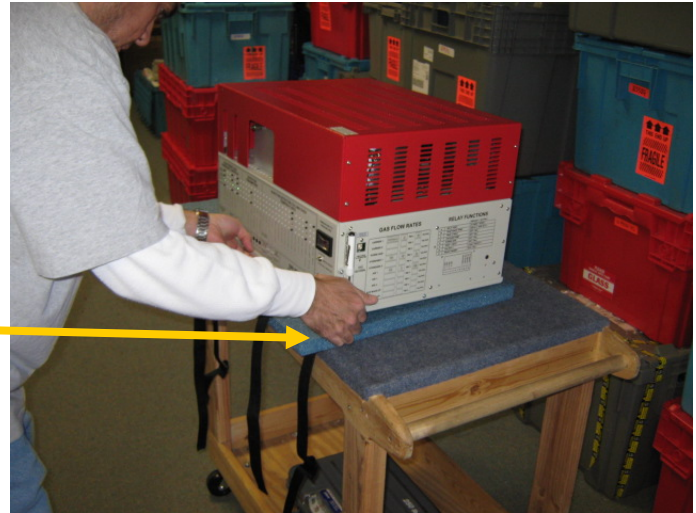

Place the black foam on top of the GC. Then put the cardboard angles on the front and back.



Tighten the two webbing straps to hold everything in place.



Moving and Packing the SRI 8610C,D or V and 310C GC for shipping

Holding the GC by the straps, lower the GC into the shipping container. The container may be grey, blue or red in color, but they are all the same size. There is room around the sides of the GC to insert accessories such as manuals, software, cables etc. There may be room for a small computer as well.

Secure the top of the container with two bolts. If the bolts are missing or lost, you can use tie-wraps or wire.

The SRI GC is now ready for shipment via FedEx, UPS or as airline baggage.

Don't forget to include your contact information inside the box in case the shipping label on the outside is lost. When sending a GC to SRI please also include a description of the problem, and a contact telephone number.

