



GC Innovations

Multiple Gas Analyzer #1

Keep your gas products in spec! Monitor gas product purity, natural gas, and ambient air quality.

Sounds expensive and complicated to operate?

Not from SRI! The SRI Multiple Gas

Analyzer #1 uses just ONE gas

sampling valve and TWO

analytical columns to perform

the same separations that

require multiple valves and

columns in other systems.

Best of all, the Multiple Gas

Analyzer #1 can achieve ppm

to 100% concentrations with a

- ★ Separates multiple gases with a single injection
- ★ Very tolerant of user adjustments and timing variations
- ★ Simpler than other multi-gas capable GC systems
- ★ Multiple gas analysis in a compact unit

19.5"W x 14.5" D x 12.5" H footprint

Heated Valve Oven

1m (3') Molecular Sieve
packed column separates
 H_2 , O_2 , N_2 , CH_4 & CO

10 Port Gas Sampling Valve
and 1mL Sample Loop

Sample In and Out
for Valve Injection

On-column
Injector

2m (6') Silica Gel packed
column separates CO_2 & C_2 - C_6

TCD Detector - universal
response, 250ppm to 100%
detection range



FID Detector - Hydrocarbon
selectivity, 5ppm detection limits
Methanizer in FID body- converts
 CO and CO_2 to Methane for FID
detection

Temperature Programmable
Column Oven

Specifically Designed for
Separation of Whole Gas Components and Natural Gas Samples

Multiple Gas Analyzer #1

Separating out the hydrocarbon components of natural gas facilitates accurate BTU quantification. This compositional analysis of a natural gas standard by an SRI Multiple Gas Analyzer #1 shows good separation up to the pentanes. Performing compositional analyses of natural gas product before and after refining helps to maximize process efficiency and profit.

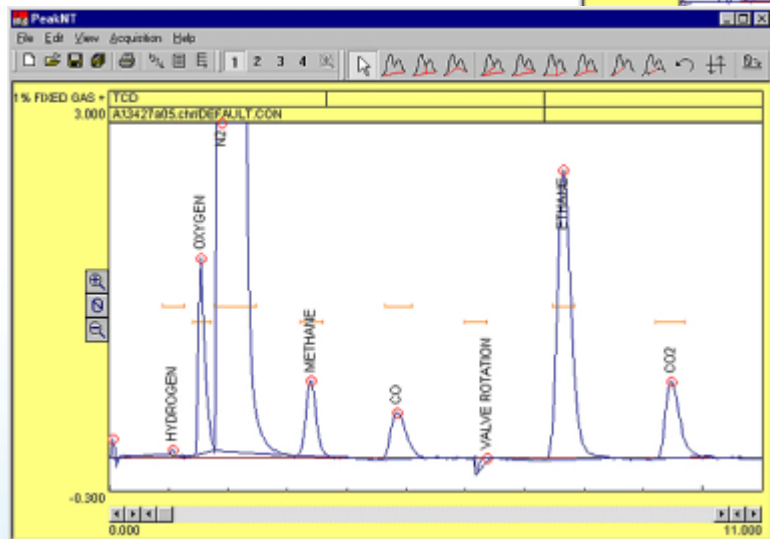
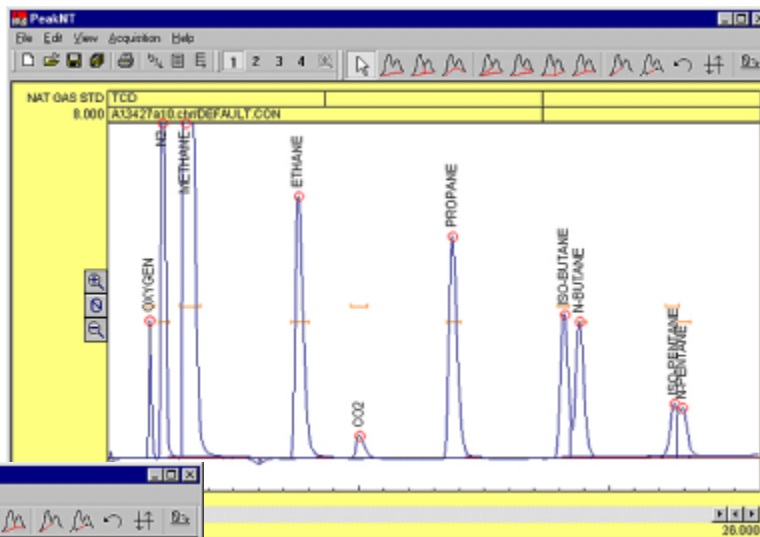


The same instrument produced this chromatogram, separating a sample mix of 1% fixed gas standard and ethane. With the built-in PeakSimple data system, the gas sampling valve was programmed to inject the sample loop contents into the carrier gas stream at 5 seconds, then rotate back at 6 minutes, after CO elution.

The basic Multiple Gas Analyzer #1 has a TCD detector only; this model provides analyses in the 250ppm to 100% range for fixed and natural gases. A second option is a TCD, Methanizer, and FID detector combination which adds 5ppm detection limits for CO, CO₂, and all hydrocarbon peaks; this model is useful for air quality monitoring and other applications. A third option is a TCD-HID detector combination, for detection limits in

the 10ppm range for all analytes...the HID even sees hydrogen! Since we build each GC from the boards up, the Multiple Gas Analyzer #1 may be further customized to suit your application needs. With the optional built-in "whisper-quiet" air compressor, the Multiple Gas Analyzer #1 can be used with the SRI H2-100 hydrogen generator to separate multiple gases anywhere, without using compressed gas cylinders!

(2022 pricing, prices subject to change, consult most recent price list.)

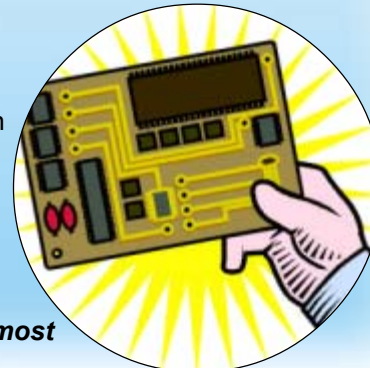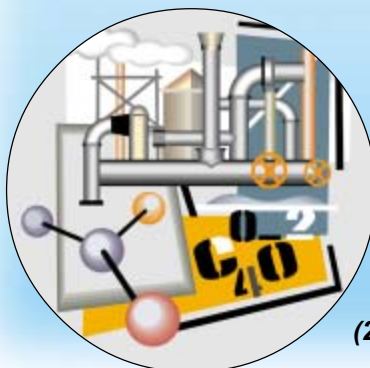

8610-0070	Multiple Gas Analyzer #1 GC with TCD detector	\$ 13,814.00
8610-0071	Multiple Gas Analyzer #1 GC with TCD, Methanizer, FID & built-in Air Compressor	\$20,029.00
8610-0072	Multiple Gas Analyzer #1 GC with TCD & HID detectors	\$20,009.00
8690-0070	Built-in Air Compressor, 120 VAC	\$ 769.00
8690-2270	Built-in Air Compressor, 220 VAC	\$ 769.00